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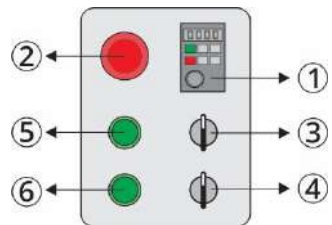
❖ **How to set up and the guideline for the laboratory mixer and permil machine:**

- Install and fix the device in an appropriate place.
- Connect the power to the 220 Voltage single phase power supply.
- Make sure the shaft and blade are not in use.
- Open the emergency power key in order to allow power to flow through the device.
- Press the lift key and lift the shaft so that the tank is completely under the mixer.
- Then lower the shaft to fit inside the tank.
- Use the lateral receivers to fully fix the tank.
- Turn the mixer switch until the shaft is ready to work.
- Use speed adjustment key to adjust the speed of the mixer shaft.

**Instructions for using the control panel:**

1. Digital Display
2. Emergency shutdown key
3. Speed adjustment key
4. Mixer key
5. Lift key
6. Down key

❖ **Service and maintenance of laboratory mixer and permil:**

- Rinse the shaft and blade after each use.
  - Grease mixer guide and main pillar.
  - Visit the power transmission belts from the electric motor to the main shaft.
  - Visit the bolts.
  - Maintenance and cleaning of the control panel and the electricity of the device.
- 
- **Please visit the above services every two months.**

# Mashinfam Company

producer of painting machines & chemical article

## Manual

Laboratory mixer and permil